

AcroSeal[®] Packaging

Your Solution for Air- and Moisture-Sensitive
Organometallic Reagents



AcroSeal Packaging Your Solution for Air- and Moisture-Sensitive Organometallic Reagents

Introduction

Organometallic reagents are widely used in organic synthesis. Typical applications include their use as strong bases, as nucleophiles, in metal exchange reactions and as reducing agents. Examples of such reagents include Grignard reagents, organolithiums, organozincs and lithium amide bases. All of these reagents react readily with moisture, oxygen or both so they must be protected under an inert atmosphere.

Historically, researchers who require organometallic reagents for synthetic purposes have chosen to freshly prepare these reagents, often *in-situ*, and use them immediately. However, safety concerns due to the highly reactive nature of these reagents, as well as a continuous push to reduce time-to-market, have led researchers to purchase commercially available organometallic reagents.

For commercially available organometallic reagents, a key concern for researchers, buyers and stockroom managers alike is how long the reagent can be stored without significant decomposition. Since these reagents are generally consumed over a period of time, the packaging system becomes even more critical.

AcroSeal packaging is specifically designed to maintain the integrity of air- and moisture-sensitive reagents while providing a convenient, easy-to-use and safe means of storing and dispensing research quantities.

Key Features of AcroSeal Packaging

Septum

AcroSeal packaging uses specially designed multi-layer septa. The AcroSeal septum used for organometallic reagents consists of a double sandwich—an inner PTFE layer that is in contact with the reagent and an outer butyl rubber layer.

The thickness of the septum is double that of typical competitor brand packaging, providing improved mechanical performance during needle puncture. The inner PTFE layer is inert to the reagent and ensures that the reagent is uncontaminated with the septum material. The thick outer butylated rubber layer allows for excellent re-sealing around needle punctures.

Repeated punctures to a septum in the same spot over time increase the risk of deformation and the likelihood that the septum will not re-seal. With competing brand packaging, it is very difficult to avoid repeated punctures in the same spot due to the very limited septum surface area. The AcroSeal septum provides more than 15 times this surface area, allowing broader spread for needle punctures, leading to better re-sealing over a longer time period.

Bottle Closure System


The AcroSeal closure system features a quadrant-style cap design that holds the septum in place, reducing the risk of bulging or deformation. A cushioned liner in the external cap presses against the septum and quadrant cap on closure, thus creating a secondary seal which prevents ingress of moisture and oxygen from the atmosphere. The tamper-evident seal, which breaks with the first opening of the cap, is used as a visible indicator of an unopened bottle.

The AcroSeal screw closure not only allows syringe withdrawal of the reagent through the septum, but provides the added convenience of permitting the top to be unscrewed and larger volumes to be poured from the bottle, which may then be closed by screwing the lid back onto the bottle.

Tamper-evident break ring at the lower end of the protective red cap. Foolproof guarantee that the customer is the first to open the bottle. The lower ring breaks off with the first opening of the red cap.

Extra thick durable rubber septum with inner PTFE resin coating. Better chemical compatibility and re-closure of any puncture hole.

Quadrant cap surface. The quadrant cap promotes the durability of the septum because punctures can be spread over the large septum surface.

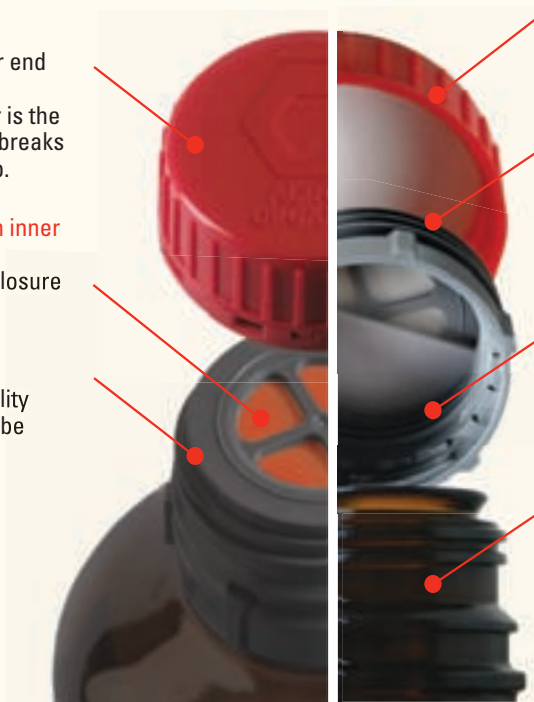

Cushioning liner in outer protective cap. Re-seals perfectly even after the septum has been punctured multiple times.

Rim design at the inner surface of the septum holding quadrant cap will hold and stretch the septum. Reduces the risk of bulging or deformation of the septum.

Three-plus windings in the septum holding cap for extra security. Large contact surface prevents infiltration of air and moisture through the screw neck.

Special rim design inside the septum holding cap fits snugly over the glass rim of the bottle neck.

No risk of accidentally unscrewing the septum holding quadrant cap.

Performance


Customers often question how well AcroSeal packaging maintains the integrity of a range of organometallics compared to the packaging utilized by other brands on the market. To address this concern, we examined the percent change in 'active' concentration of a range of organometallics stored in AcroSeal packaging and in competing brand packaging.

Test Results and Discussion

An eight-week study of 'active' concentration of four organometallic reagents was conducted to determine the effectiveness of the AcroSeal packaging system and a competitor's packaging in maintaining the integrity of the organometallic reagent under normal laboratory usage.


n-Butyllithium, 1.6M in hexanes

Over an eight-week period, the change in concentration of *n*-Butyllithium in hexanes in the AcroSeal packaging was negligible within the experimental error. Over the same time frame, *n*-Butyllithium stored in the competitor's packaging showed more than a 15% drop in 'active' concentration.

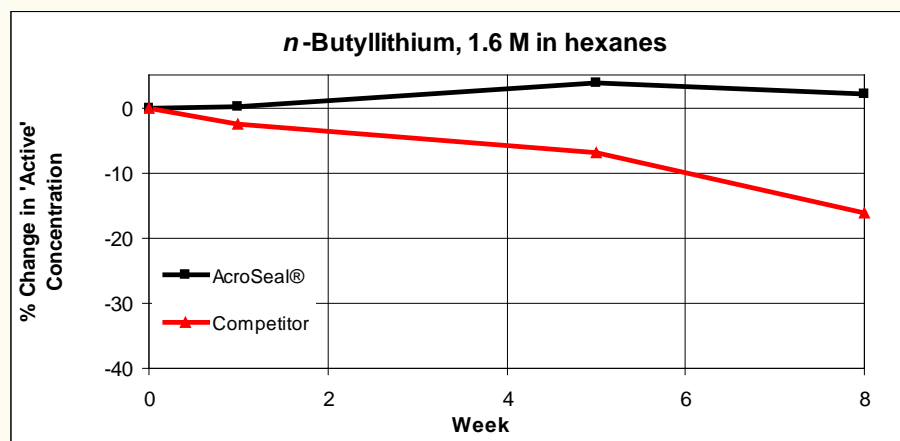


Phenylmagnesium chloride, 2M in tetrahydrofuran

Up until week 5, the 'active' concentration of Phenylmagnesium chloride, 25 wt% (2M) in tetrahydrofuran, stored in AcroSeal packaging remained constant (within experimental error), with a less than 5% drop by week 8. The 'active' concentration in the competitor product dropped rapidly by around 10% after the first week and lost more than 15% activity by week 8.

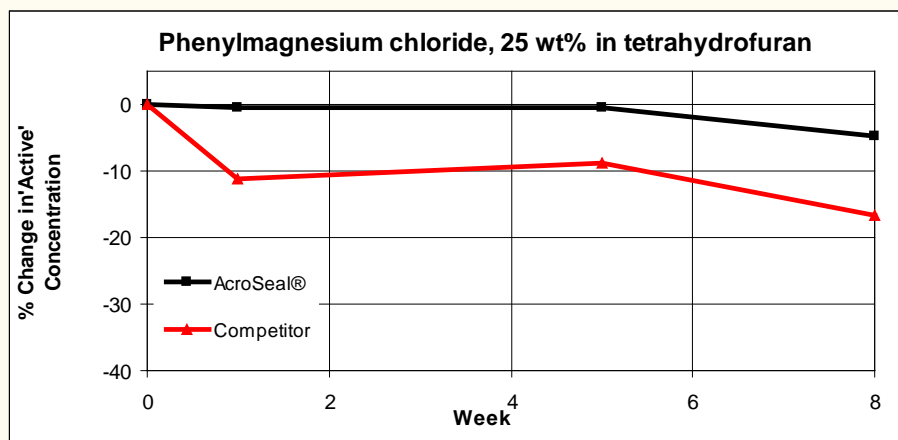


Ethylmagnesium bromide, 1M in tetrahydrofuran

For Ethylmagnesium bromide, 1M in tetrahydrofuran, the rate of loss of active organometallic was found to be twice as fast with the competitor's packaging than with AcroSeal. After eight weeks, the active level decreased by 50% in the competitor's packaging.

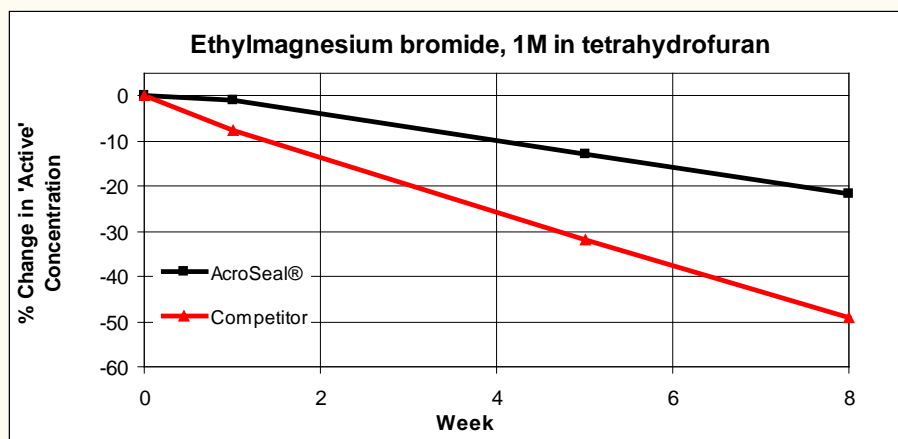

LDA, 2M in tetrahydrofuran/*n*-heptane/ethylbenzene

Lithium diisopropylamide, 2M in tetrahydrofuran/*n*-heptane/ethylbenzene was relatively stable in both AcroSeal and competitor packaging over a five-week period. After eight weeks the 'active' concentration dropped by 4% in the AcroSeal packaging and 11% in the competitor packaging.

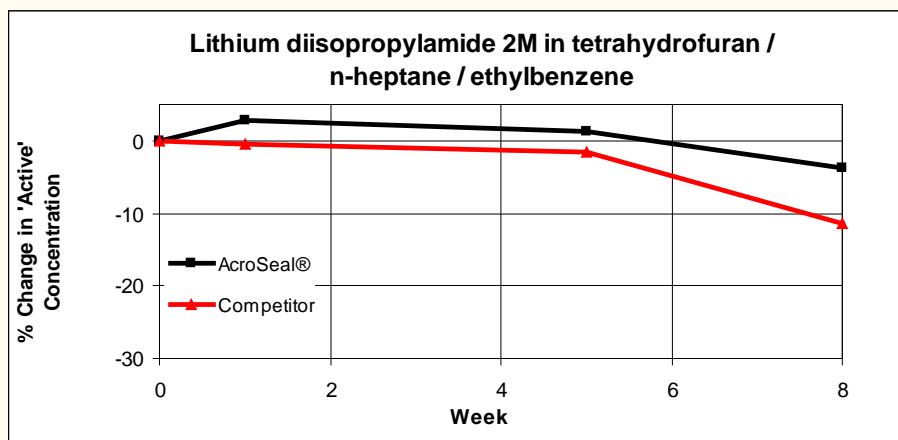



Conclusions

This study has shown that AcroSeal packaging is suitable for a wide variety of organometallics dissolved in a range of solvents:

- *n*-Butyllithium, 1.6M in hexanes
- Phenylmagnesium chloride, 25 wt% (2M) in tetrahydrofuran
- Ethylmagnesium bromide, 1M in tetrahydrofuran
- Lithium diisopropylamide, 2M in tetrahydrofuran/*n*-heptane/ethylbenzene

In all cases, the AcroSeal packaging maintained the integrity of these organometallic reagents handled under normal laboratory conditions over several weeks better than the competitor packaging.

For a complete list of reagents available in AcroSeal packaging, please visit our website at www.acros.com.

Experimental Section

Introduction

Solutions of organometallic reagents in inert solvents inevitably contain some metal oxides, hydroxides, alkoxides (or aroxides) and even carbonates, which result from reaction with oxygen, moisture and atmospheric carbon dioxide. If a septum fails to re-seal properly over time, the concentration of the 'active' organometallic species decreases, while the concentration of the inorganic basic by-products increases. In these studies, 'active' concentration refers to the concentration of the organometallic species.

The 'active' concentration of *n*-butyllithium in hexanes was measured using the Gilman double titration method¹, while the Watson-Eastham method was used for the other reagents².

Samples were taken over eight weeks, in duplicate, from 100mL bottles. The initial measurement is designated as week zero. Two bottles of each reagent with identical batch numbers were analyzed and the four results were averaged. Using 21-gauge needles, each bottle received five new punctures per septum, per week of the study. At the end of each test, the caps were firmly secured by hand. The organometallic reagents were stored and analyzed at ambient temperature.

Packaging performance was compared by examining the percent change in 'active' concentration over time. The initial 'active' concentration was subtracted from successive testing values to normalize the concentration over time. Division by the initial value and multiplication by 100 gives the percent change in 'active' concentration.

General Handling

50 mL round-bottom flasks, magnetic stirrer bars and B19 rubber septa were dried overnight in an oven at 110°C. They were removed and the flask, containing a stirrer bar, was sealed with the B19 septum. The flask was evacuated under vacuum and filled three times with nitrogen gas via a manifold.

Disposable syringes and needles used during the procedures were flushed with nitrogen several times immediately prior to use.

During dispensing of the reagent, a nitrogen line from the manifold was attached to each bottle via a syringe needle to replace the volume of reagent drawn up in the syringe.

Double Gilman Titration Method

Determination of the total base content:

A nitrogen-flushed syringe was filled with approximately 5mL of the *n*-butyllithium in hexane solution and weighed (W1). The reagent was injected into a round-bottom flask containing anhydrous hexanes (20mL) under a nitrogen atmosphere. The empty syringe was then weighed (W2). Water (20mL, HPLC grade) was added and the mixture was stirred for 10 minutes. Phenolphthalein indicator (2 drops) was added and the mixture was titrated against 1.0M hydrochloric acid (HCl) until complete disappearance of the pink color.

$$\text{Calculation: Percent total base} = \frac{(V \times C)}{[(W1 - W2)/D]}$$

Where V = titer of HCl added (mL)

C = concentration of HCl (M)

W1 = weight of syringe containing *n*-butyllithium (grams)

W2 = weight of syringe after *n*-Butyllithium addition (grams)

D = density of the *n*-Butyllithium solution (0.68g/mL 25°C)

Determination of the free base content:

A nitrogen-flushed syringe was filled with approximately 5mL of the *n*-butyllithium in hexane solution, and weighed (W1). The reagent was injected into a round-bottom flask containing anhydrous hexanes (20mL) under a nitrogen atmosphere, followed by the addition of anhydrous diethyl ether (0.2mL). The empty syringe was then weighed (W2). Freshly distilled benzyl chloride (5mL) was added dropwise and the mixture was stirred for 15 minutes to ensure complete reaction with the 'active' organometallic. Water (25mL, HPLC grade) was added and the mixture was stirred for 10 minutes. Phenolphthalein indicator (2 drops) was added and the mixture was titrated against 0.1M HCl until complete disappearance of the pink color.

$$\text{Calculation: Percent free base} = \frac{(V \times C)}{[(W1 - W2)/D]}$$

Calculation of the 'Active' base content:

Percent 'active' base = Total base – Free base

References

¹Gilman, H.; Haubein, A. H.; *J. Am. Chem. Soc.*, 1944, 66, 1515

²Watson, S. C.; Eastham, J. F.; *J. Organometallic Chem.*, 1967, 9, 165



Watson-Eastham Method

A nitrogen-flushed syringe was filled with approximately 5mL of the organometallic reagent solution and weighed (W1). This was injected into a round-bottom flask containing anhydrous toluene (20mL) and approximately 8mg of the appropriate indicator under a nitrogen atmosphere. The empty syringe was then weighed (W2).

1,10-Phenanthroline was used as an indicator for the Grignard solutions. The mixture was titrated against freshly prepared and standardized approximately 0.1M *sec*-butanol solution in anhydrous toluene until disappearance of the pink color.

2,2'-Biquinoline was used as an indicator for the lithium diisopropylamide solution. The mixture was titrated against freshly prepared and standardized approximately 0.1M *sec*-butanol solution in anhydrous toluene until disappearance of the red color.

Calculation of the 'active' base content:

$$\text{Calculation: \% 'Active' base} = \frac{(V \times C)}{[(W1 - W2)/D]}$$

Where V = titer of *sec*-butanol added (mL)

C = concentration of *sec*-butanol (M)

W1 = weight of syringe containing the organometallic reagent (grams)

W2 = weight of syringe after organometallic reagent addition (grams)

D = density of the organometallic solution (g/mL, 25°C)

Densities of the organometallic reagents:

Phenylmagnesium chloride, 25 wt% (2M) in tetrahydrofuran (d=1.042g/mL, 25°C)

Ethylmagnesium bromide, 1M in tetrahydrofuran (d=1.02g/mL, 25°C)

Lithium diisopropylamide, 2M in tetrahydrofuran/*n*-heptane/ethylbenzene (d=0.812 g/mL, 25°C)

GLOBAL LOCATIONS
AMERICAS
Canada

Fisher Scientific Canada
112 Colonnade Road
Ottawa, Ontario
Post Code: K2E 7L6
Toll-Free Number: 800-234-7437
Fax: 800-483-2996
www.fishersci.ca

Latin America

Fisher Scientific Global Export, Latin America
2970 Johns Creek Court
Suite 500
Sawnee, GA
Post Code: 30024
Toll-Free Number: 770-871-4725
Fax: 770-871-4726
www.fishersci.com

United States

Fisher HealthCare
9999 Veterans Memorial Drive
Houston, TX
Post Code: 77038
Toll-Free Number: 800-640-0640
Fax: 800-290-0290
www.fisherhealthcare.com
Fisher Scientific
2600 Park Lane Drive
Pittsburgh, PA
Post Code: 15275
Toll-Free Number: 800-756-7000
Fax: 800-926-1166
www.fishersci.com

ASIA
China

Fisher Scientific China
Toll-Free Number: 400 881 5117
sales.china@thermofisher.com
www.fishersci.com.cn

Shanghai Corporate Office
6/F Long Life Mansion
No. 1566 Yan An West Rd.
Shanghai, China
Post Code: 200052
Tel: (8621) 5258 1100
Fax: (8621) 5258 0119

Beijing Office
Units 707-715, 7th Floor
Tower West, Yonghe Plaza
No. 28 Andingmen East Street
Beijing, China
Post Code: 100007
Tel: (8610) 8419 3588
Fax: (8610) 8419 3580

Guangzhou Office
Room 2405-2406, JianLiBao Mansion
No. 410-412 Middle DongFeng Rd.
Guangzhou, China
Post Code: 510030
Tel: (8620) 8214 5288
Fax: (8620) 3877 1941

India

Fisher Scientific India
403-404, B-Wing, Delphi
Hiranandani Business Park
Powai, Mumbai - 400 076
Tel: 1 800 208 7001
Fax: +91 22 6680 3001/3002
qf.customercare@thermofisher.com
http://www.fishersci.in

Japan

Fisher Scientific Japan
Thermo Fisher Scientific K.K.
C-2F, 3-9 Moriya-cho
Kanawaga-ku, Yokohama
221-0022 Japan
Tel: 81 45 450 6310
Fax: 81 45 450 6318
support@fishersci.jp
www.fishersci.co.jp

Korea

Fisher Scientific Korea
Sambu Bldg. 13F
676 Yeoksam-dong,
Kangnam-Gu
Seoul 125-879, Korea
Dir: (02) 527-0300
Customer Service Center: (02) 527-0300
Fax: (02) 527-0311
sales@fishersci.co.kr
www.acros.co.kr

Malaysia

Fisher Scientific Malaysia Sdn Bhd
No. 3 Jalar Sepadu 25/123
Taman Perindustrian Axis Sekayen 25
40400 Shah Alam
Selangor Darul Ehsan Malaysia
Technical Service Hotline: 1-300-88-7888
Tel: (603) 5'218888
Fax: (603) 51218899
enquiry.my@thermofisher.com
www.fishersci.com.my

Singapore

Fisher Scientific Singapore
Fisher Scientific Pte Ltd.
8 Pandan Crescent
LLA, #05-05 UE Tech Park
Singapore 128464
Tel: (65) 6873 6006
Fax: (65) 6873 5005
enquiry.sg@thermofisher.com
www.fishersci.com.sg

EUROPE
Austria

Fisher Scientific (Austria) GmbH
Rudolf von Alt-Platz 1
A-1030 Wien
Phone: 0800 20 88 40
Fax: 0800 20 66 90
info.austria@thermofisher.com
www.at.fishersci.com

Belgium

Fisher Scientific
BP 567
B-7500 Tournai 1
Tel: 056 260 260
Fax: 056 260 270
be.fisher@thermofisher.com
www.be.fishersci.com

Czech Republic

Fisher Scientific, spol. s r.o.
Kosmonautů 234
Pardubice
CZ-530 09
Tel: 466 798 230
Fax: 466 435 008
info.cz@thermofisher.com
www.thermofisher.cz

Denmark

Fisher Scientific Biotech Line A/S
Industrivej 3
Postboks 60
DK-2650 Slangerup
Tel: +45 70 27 99 20
Fax: +45 70 27 99 29
kundeservice@thermofisher.com
www.fishersci.dk

Finland

Fisher Scientific Oy
Ratasitie 2
FI-01620 Vantaa
Tel: +358 902 76 237
Fax: +358 902 76 235
fisher.fi@thermofisher.com
www.fishersci.fi

France

Fisher Scientific
Parc d'Innovation BP 50111
67403 Illkirch Cedex
Tel: 03 88 67 53 20
Fax: 03 88 67 11 68
fr.commande@thermofisher.com
www.fr.fishersci.com

Germany

Fisher Scientific GmbH
Im Heiligen Feld 17
D-58239 Schwerte
Phone: 0800 3 47 43 70
Fax: 0800 3 47 43 71
info.germany@thermofisher.com
www.de.fishersci.com

Ireland

Fisher Scientific Ireland
Suite 3 Plaza 212
Blanchardstown Corporate
Park 2
Ballycoolin
Dublin 15
Tel: +353 01 885 5054
Fax: +353 01 899 1855
ie.sales@thermofisher.com
www.ie.fishersci.com

Italy

Fisher Scientific
Tel: 02 953 26 258
Fax: 02 953 27 274
it.fisher@thermofisher.com
www.it.fishersci.com

The Netherlands

Fisher Scientific
Postbus 4
1120 AA Landsmeer
Tel: 020 487 70 00
Fax: 020 487 70 70
nl.info@thermofisher.com
www.nl.fishersci.com

Norway

Fisher Scientific
Frystaveien 33E
0884 Oslo
Tel: +47 22 95 53 58
Fax: +47 22 95 59 40
fisher.no@thermofisher.com
www.fishersci.no

Portugal

Fisher Scientific
Rua Pedro Álvares Cabral, nº24, 2ºD
Edifício Euro - Infanteado
2670-391 Loures
Telephone: +351 21 425 33 50/4
Fax: +351 21 425 33 51
pt.fisher@thermofisher.com
www.pt.fishersci.com

Spain

Fisher Scientific
C/ Luis I, 9
28031 MADRID
Tno: 91 380 87 10
Fax: 91 380 85 02
es.fisher@thermofisher.com
www.es.fishersci.com

Sweden

Fisher Scientific
Box 3193
400 94 Göteborg
Tel: +46 31 - 68 94 30
Fax: +46 31 - 68 07 17
gt.info@thermofisher.com
www.fishersci.se

Switzerland

Fisher Scientific
Wilstrasse 57 - Postfach 1000
5610 Wohlen
Tel: 056 618 41 11
Fax: 056 618 41 41
info.ch@thermofisher.com
www.ch.fishersci.com

United Kingdom

Fisher Scientific UK
Bishop Meadow Road
Loughborough
Leicestershire LE11 5RG
Tel: +44 (0)1509 231186
Fax: +44 (0)1509 231803
info@fisher.co.uk
www.fisher.co.uk

Rest of Europe

Thermo Fisher Scientific
Geel West Zeele 2
Janszaan Pharmaceuticaan 2a
2440 Geel - Belgium
Tel: +32 34 57 52 11
Fax: +32 14 59 26 10
www.acros.com

MIDDLE EAST AND AFRICA
Fisher Scientific Global Export,

Latin America
2970 Johns Creek Court
Suite 500
Sawnee, GA
Post Code: 30024
Toll-Free Number: 770-871-4725
Fax: 770-871-4726
www.fishersci.com

OCEANIA
Australia

Thermo Fisher Scientific Australia Pty Ltd
5 Caribbean Drive
Scoresby, VIC 3179
Tel: +1300 735 292
Fax: +61 3 9763 1169
AUinfo@thermofisher.com
www.thermofisher.com.au

New Zealand

Thermo Fisher Scientific New Zealand Ltd 244
Bush Road,
Albany, Auckland 0632
Tel: 0800 833 966
Fax: +64 9 580 6788
info.nz@thermofisher.com
www.thermofisher.co.nz

©2010 Thermo Fisher Scientific Inc.
All rights reserved. All brands and trademarks
are part of Thermo Fisher Scientific Inc. and
its subsidiaries.