

Thermo Scientific
ALPS 50V
User's Guide



Important:

IT IS ESSENTIAL THAT THE USER OF THIS MANUAL IS AWARE OF THE POTENTIAL HAZARDS ASSOCIATED WITH THE UNIT AND ITS ACCESSORIES.

ALL OPERATORS SHOULD BE FAMILIAR WITH THE SAFETY PRECAUTIONS AND WARNINGS GIVEN IN THIS SECTION PRIOR TO ATTEMPTING TO OPERATE THE UNIT.

IF THE UNIT IS USED IN A MANNER NOT SPECIFIED BY THE MANUFACTURER, THE PROTECTION PROVIDED BY THE EQUIPMENT MAY BE IMPAIRED.


The following symbols are used in this manual and on the instrument:




CAUTION/WARNING



HEATED SURFACE

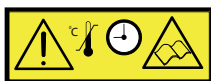

Refer to manual to set correct time and temperature

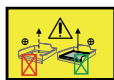

Ensure correct adapter plate is fitted before operating

The information in this user guide including any apparatus, Methods, Techniques and Concepts described herein, are the property of Thermo Fisher Scientific or its Licensees and may not be copied, disclosed or used for any purpose not expressly authorized by the owners thereof.

Since Thermo Fisher Scientific constantly strives to improve all of its products, we reserve the right to modify equipment and any user instruction manual without prior notice. No part of this instruction manual may be reproduced in any form without the prior consent of Thermo Fisher Scientific.

Thermo Fisher Scientific
22 Friars Drive
Hudson NH, 03051

© 2007 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries.



WARNING: If the unit is used in a manner not specified by the manufacturer, the protection provided by the equipment may be impaired.

- The unit is intended for indoor laboratory use only, at an altitude of less than 2200m above sea level, within a temperature range of 18°C to 30°C and a relative humidity range of 20% to 80% non-condensing. If the instrument is stored outside these ranges, it should be left to stand until it equilibrates to within the above limits.
- Ensure the voltage selection switch is set to the correct voltage for the required region (See Fig 1)
- Ensure the correct fuse for the required voltage setting is fitted (see Specifications section).
- Do not operate the unit outside the rated power supply range specified.
- There are no user accessible or serviceable parts inside the unit. Do not remove or open the unit's casework.
- Before using any cleaning or decontamination method except those recommended by the manufacturer, users should check with the manufacturer that the proposed method will not damage the equipment.
- Ensure the unit is only connected to an earthed supply.
- Ensure adapter plate is fitted before use.
- Once timer starts, no additional handle pressure is required.
- Adjust time and temperature only for optimum seal.



WARNING: The hot plate can reach temperatures up to 200°C. Care must be taken not to touch it or serious burns may occur.

- The unit will remain hot for a considerable time after it has been turned off and must be allowed to cool down to an acceptable level before cleaning.
- Foil sealed plates may remain hot for a number of seconds after sealing and should be handled with care

Contents

General Description of Apparatus	Page 5
Intended Function of Apparatus.....	Page 5
Installation.....	Page 5
Types of Consumables.....	Page 6
Operating Instructions	Page 6
Operating Instructions	Page 7
Operating Instructions:	Page 8
Cleaning.....	Page 8
Specifications	Page 8
Over Temperature Safety Cutout	Page 9
Switching off the Audible Warning	Page 9
Lifting and moving the instrument.....	Page 9
Warranty Returns	Page 9
Declaration of Conformity	Page 10

General Description of Apparatus

The Thermo Scientific ALPS 50V provides a safe, controlled method of sealing sample plates made from various plastics. The sealer is adjustable for both sealing time and temperature and gives a visual countdown display during the sealing process. The ALPS 50V can seal deep well plates (Bioblocks) with various thermal sealing plastic or foil films and may be adapted to seal Microplates and PCR plates by using a suitable plate carrier.

Intended Function of Apparatus

The ALPS 50V is intended to provide a platform with which to load and thermally seal biological plates to prevent sample loss at high temperature or prolonged periods of storage. Various types of plate may be loaded onto a plate carrier and covered by a thermal sealing medium, which may be foil or clear plastic. The operator may pre-set a precise temperature and seal time for the sealing operation. Sealing is initiated by pulling down a handle which raises the sample plate into contact with the hot plate. The plate sealer detects when the sample plate is in the correct position and starts a countdown timer. The countdown is displayed on a three digit LED display. When the display reads zero, the operator lifts the handle back to the upright position, which moves the sample plate away from the hot plate. The plate may then be manually removed for further processing.


Installation

The variable temperature sealer should be removed from its packaging and all internal packaging removed. The instrument should be placed on a level surface, away from direct sunlight and drafts ensuring access to the power switch on the rear of the unit. Ensure that the vents on the back panel are not obstructed.

Mains inlet selection:

The mains voltage selector switch should be checked to ensure that the desired voltage has been set (Fig 1) and the correct value of fuse has been fitted (see FIG 2 and the specification section).

Fig 1

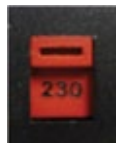

Voltage Selector Switch Set to 230V



Voltage Selector Switch Set to 115V

Fig 2



Fuse Position

Fig 3



Main Power Switch

Types of Consumables

The ALPS 50V is designed to be used with the following consumables:

Plate Types: (Polypropylene, polyethylene or polystyrene)

- Deep well plates without a carrier.
- PCR (Polymerase chain reaction) plates with carrier AB-0563/1000 and AB-0724.
- Microplate with carrier AB-0827.

Sealing film types:

- AB-0559 Thermo Seal
- AB-0757 Easy Pierce
- AB-1720 Easy Piece 20µm
- AB-0745 Easy Peel
- AB-1797 Clear Seal
- AB-0812 Clear Seal Diamond
- AB-0685 Clear Seal Strong

Product Includes

- 96 well PCR plate adapter (AB-0563/1000)
- BioBlock Adaptor Plate
- US Power Cord
- EU Power Cord
- UK Power Cord
- 3.15A Fuse
- User Guide Manual: For translations in French, Italian, Danish, Spanish and Portuguese please go to: XXXXXXXXXXXXXXXXXXXX

Package Note

M9800119: Registration Form ALPS 50V

Operating Instructions

If the instrument has been removed from storage, it should be left to reach ambient room temperature. For correct operating conditions, please see specification section. Please refer to (Fig 5) for location of user interface components and (Fig 3) for the location of the power switch. When the instrument is first turned on, The LED display shows the current settings for either sealing temperature or seal time and the heater on/off status retained from the last seal. Pressing the mode switch toggles the display between seal time and temperature setting. A new temperature setting between 125°C and 200°C may be set by pressing the up and down arrow keys. The “C” mode LED lights to indicate that the adjustable field is temperature. Seal time is adjustable via the up and down keys in 0.5 Second increments between 1 – 9 seconds. If the up or down arrow keys are held for more than 1 second, the key will auto-repeat. When the desired settings have been set the heat on/off button may be pressed to begin heating the sealing plate to the set temperature. The heat LED flashes at a slow rate to indicate that the heater is on but not yet up to temperature. The heater should heat up to the set point in about ten minutes. When the heater is within two degrees of the set point, the heat LED will be continuously on and the sealer is ready to begin sealing plates.

Operating Instructions

A Typical operating sequence is:

1. Switch instrument on and set required sealing time and heater set-point temperature for the desired sealing process, via the keypad and LED display.
2. Enable the heater by pressing the Heat on/off button.
3. Allow sufficient time for the instrument to reach desired temperature. The Heat on/off LED will flash whilst the sealer is coming up to temperature and will remain on permanently when the set point ($\pm 2^{\circ}\text{C}$) is reached.
4. Load sample-plate onto plate carrier and add the sealing film. Care should be taken not to touch the heating surface whilst loading the sample plate.
5. Grasp the handle with one or two hands as appropriate (see FIG 4a). When appropriate i.e. if one hand is used, ensure free hand is away from heat source and moving parts.
6. Lower handle to thermally compress sealing film onto sample plate (see FIG 4b).
7. When correct clamping pressure is achieved an audible warning will sound (if audible sounder fitted) and the “timer” will count down to zero. Care should be taken not to apply more pressure to the handle than is necessary to operate the micro switch.
8. On zero seconds, the display will flash and an audible warning will sound. Release the heater plate from the sealed sample plate by raising the handle to its upper position and review the seal that has been made. Failure to raise the handle could ruin the sample plate.

Note: The heater does not turn off when the buzzer stops sounding

9. Remove sealed plate for subsequent processing.

The seal temperature and time may vary with each type of plate. It is recommended that the temperature and time settings for the sealer are reviewed empirically to determine optimum seal integrity. The final determination of whether a seal has been made or not is dependent on the users processes

Operating Instructions:

Fig 4a



Fig 4b

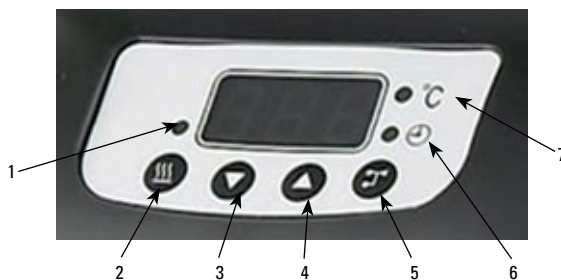


Fig 4c



Fig 5

1. Heat On/Off LED
2. Heat On/Off Button
3. Down Button
4. Up Button
5. Mode Button
6. Set Seal Time Mode LED
7. Set Temperature Mode LED



Cleaning

The instrument is to be cleaned only by wiping with Virkon solution at the manufacturer's guideline concentration. No parts are to be autoclaved. Failure to comply with the manufacturers requirements may cause damage to the function and operation of the instrument.

Specifications

Equipment Reference:
ALPS 50V™

Model No:
953-A

Specification:

Temperature Set Point Range	125°C to 200°C in 1°C increments.
Time Set Range	1 to 9 seconds in 0.5 second increments.
Operating Humidity Range	20% to 80% non-condensing.
Operating Temperature	18°C to 30°C.
Size	W 220 x D 321 x H 425 mm
Weight	7.2 kg
Range	100 - 130V AC 50/60Hz or 220 – 240V AC 50/60Hz
Power Supply	350 W Max
Power Rating	100 - 130V - T3.15A (IEC 127) or
Fuse	220 - 240V - T1.6A (IEC 127)
Mains Input Connector	IEC 320
Applicable Standards	FCC Class A, CFR 47 Part 15 BS EN 61326:1998

Over Temperature Safety Cutout

The instrument is protected against a fault condition leading to the sealing plate heating up to a temperature exceeding its maximum set point. Protection is provided by a single use thermal cut out, operating in the range 235°C to 240°C. In the unlikely event that this cut out should operate, it would necessitate the sealer being returned to the manufacturer for repair.

Switching off the Audible Warning

Some units are fitted with a sounder which give an audible warning at the end of the seal time. The sounder is on by default but may be programmed to be on or off. To turn the Audible warning off, place the sealer in Temperature set mode (Set Temperature Mode LED on) and simultaneously press the Mode button and the Heat On/Off buttons together (See Fig. The display will show "01" to indicate that the audible warning is ON and both Set Temperature Mode and Set Seal Time Mode LEDs will be off. Pressing the Down button turns the audible warning OFF. The display should now show "00". Pressing the Mode button will save this configuration so that the audible warning is always off and the display will return to the Set Temperature Mode. To turn the audible warning back on repeat this process but instead press the Up button to display "01".

Lifting and moving the instrument



WARNING: Before attempting to lift the instrument after use it should be left for a period of 30 minutes to allow the dissipation of residual heat in the heater pad, care should be taken not to touch any of the heated surfaces even after the cooling period. The instrument should be lifted by holding the operating handle. Care should be taken when lifting the instrument due to its weight. Refer to current Health and Safety procedures.



← Operating handle

Warranty Returns

Thermo Fisher Scientific
22 Friars Drive
Hudson NH, 03051
(800) 345-0206
www.thermo.com
The world leader in serving science

Returns will not be accepted without prior authorization. Prior to the return of equipment please contact Technical Service at 1-800-345-0206; all equipment that may be contaminated with biohazardous materials, potentially biohazardous materials or radioactivity, must be decontaminated. A form will be sent to you and it must be returned to Matrix Technologies part of Thermo Fisher Scientific with your instrument or by faxing it to (603) 595-0106.

Declaration of Conformity

The ALPS 50V has been designed in accordance with, and satisfies the requirements of, article 11 of the Low Voltage Directive 73/23/EEC as realigned by 93/68/EEC on the harmonization of the laws of the Member States relating to electrical equipment designed for use within certain voltage limits, to the essential requirements of BS EN 61010-1:2001.

The ALPS 50V has been type tested by EMC Projects Limited (a UKAS and CAA approved test facility and UK appointed Notified Body), and issued a Certificate of

Compliance No. 6091/06 to the following EMC standard:

BS EN 61326:1998, Electrical equipment for measurement, control and laboratory use. EMC requirements

Satisfying the EMC Directive(s) 89/336/EEC and 92/31/EEC as realigned by 93/69/EEC.

This instrument has been tested and meets the requirements of the following CSA classes;

CLASS 8721 04 – LABORATORY ELECTRICAL EQUIPMENT

CLASS 8721 84 - ELECTRICAL EQUIPMENT FOR LABORATORY USE (Certified to U.S. Standards) and as such bears the CSA logo

Brian Archbold
Director Operations

Thermo Fisher Scientific
Liquid Handling Consumables
22 Friars Drive
Hudson, NH 03051 USA

Thermo Fisher Scientific
22 Friars Drive
Hudson NH, 03051
(800) 345-0206

The world leader in serving science

© 2010 Thermo Fisher Scientific Inc. All rights reserved.
All other trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries.

www.thermoscientific.com

North America: Tel: 800 345 0206 | email: matrix.info@thermofisher.com
Europe: Tel: +44 (0) 161 486 2110 | email: matrix.eu.info@thermofisher.com
Asia: China: Tel: +86-21-6865 4588, **Japan:** Tel: +81-45-453-9220, **India:** +91-22-6716-2200,
Hong Kong & Other Asian Countries: Tel: +852-2885-4613, **Australia:** Tel: 1-300-735-292,
New Zealand: Tel: 0800-933-966 | email: matrix.ap.info@thermofisher.com

Rev A

Thermo
SCIENTIFIC